


# BEFORE YOU BEGIN

Please remember that the 2100 Retriever is an autoclave and should be handled appropriately.

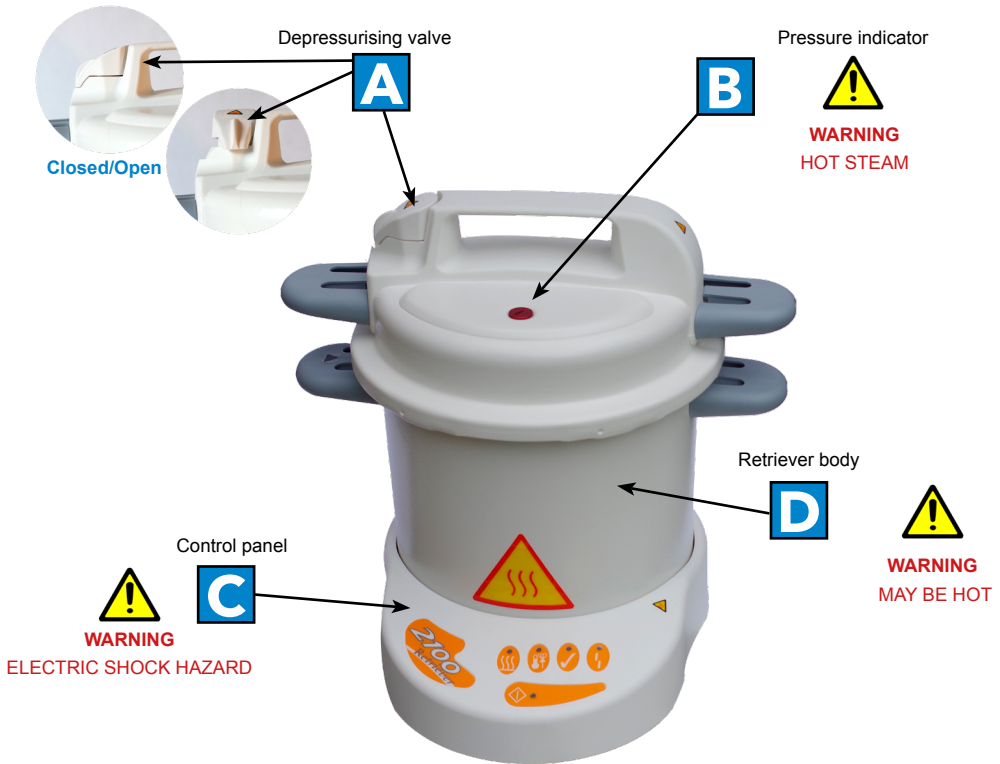

## DO NOT DO THE FOLLOWING:

- Touch the unit whilst in operation - it gets hot.
- Attempt to remove the lid during operation.
- Lubricate the gasket.
- Add any chemicals to the water poured into the unit.
- Attempt to use volatile substances and toxic materials in buffers, or inappropriate loads such as easy melting plastics.
- Place the unit on heat sensitive surfaces (polished wood or glass).
- Open the depressurisation valve (A) during the cycle.
- Press the start button once the cycle has been started, as this will reset the cycle.
- Leave the depressurisation valve (A) in the "Open" position when placing the lid upside down on a work surface.
- Immerse the unit or electrical cord in water when cleaning.
- Use in areas of risk associated with flammable materials or gasses.
- Attempt to change fuses until the unit has been unplugged from the mains. Only qualified persons should change fuses.
- Reach over the unit when removing the cover, to do so may cause burns from rising heat and steam.